# Choose your language

# PP-Inspector

# **Anleitung**

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PC-Software for the PP-Serie Power-Potentiostats


### 1. Installation

During the installation of the THALES- Software, a folder 'C:\THALES\PPINSPECTOR' is created containing the files required to run a power potentiostat of the PP series (PP- device).

To run a PP- device, start the program *PP2xx.exe*. Executing this program for the first time, the following message is displayed on the screen:

Click on 'Yes' for a first time installation. Now, for the creation of an initialization file, the software analyses your system (requires about 5 seconds) and the following screen will be displayed: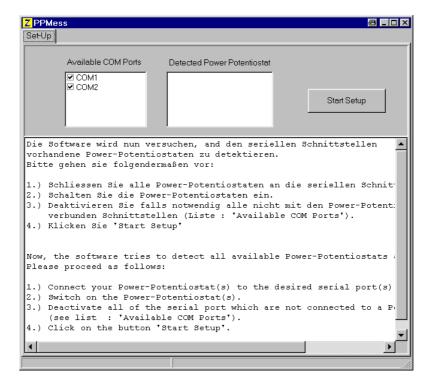

In the table **Available Comports** all available COM ports of the system are listed. The table **Detected Power Potentiostats** is still empty. Clicking on the button on the right side (**Start Setup**) will start the setup procedure.

Note: The software is able to handle more than one power potentiostat at the same time. Provided you want to add another potentiostat later, just delete the file **PP200.ini** which is created during the initialization procedure and run the program.


## 2. Settings

In this section of the software, the static parameters of the PP- device are set:

- 6. **Select Power Potentiostat :** Provided you have more than one PP- devices connected to your PC, you can set one of them to the active PP- device. Note: *type* (PP200, PP210 or PP240) as well as the involved COM- port are displayed in the header of the window too
- 7. **Shunt**: To select a shunt resistor
- 8. **Bandwith**: A property of special interest for impedance measurements. Keep the value of 10 KHz.
- 9. **Full Scale**: Here you set the maximum measurable potential. With respect of best resolution, try to keep the default setting  $\pm 4$  V if possible.
- 10. **Mode:** changes between potentiostatic and galvanostatic mode of operation.

On the left hand side the actual measured values of potential and current are displayed. Below, you can insert the value to be set. You can use different notations to set this value, i. e. Entries of 0.1 or 0,1 or 1e-1 or 100m will set a value of 100 mV in potentiostatic and 100 mA in galvanostatic mode of operation.

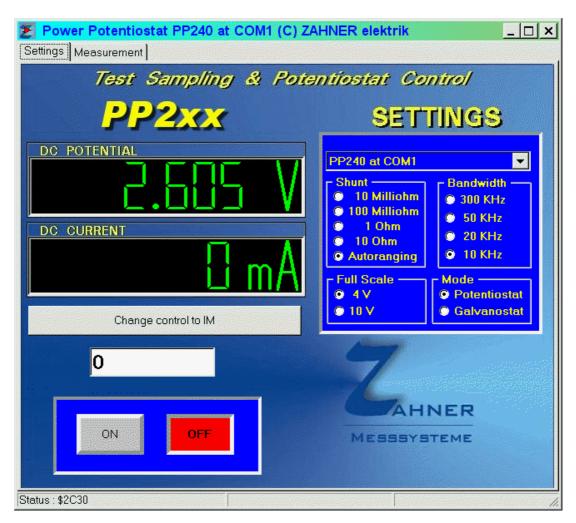
Clicking on the button **Change control to IM** / **Change control to PC**, you can toggle between these two devices to control the Power Potentiostat.

### 3. Measurement

In this menu you can set/select static as well as dynamic parameters for measurements depending on the radio button **Mode of measurement**.

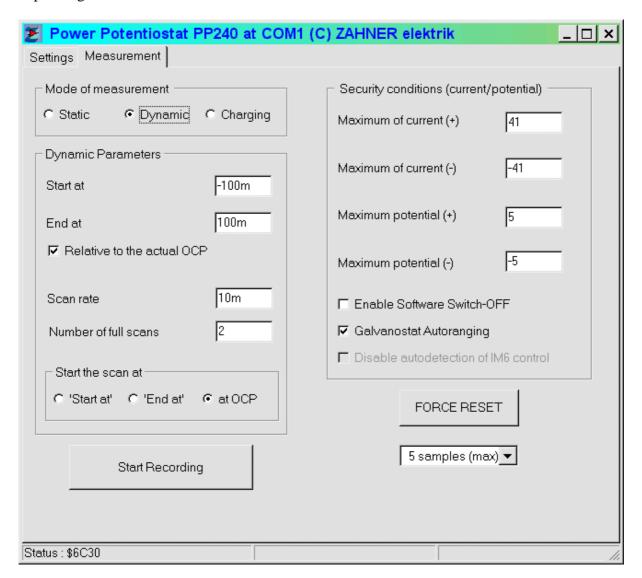Independent of static or dynamic mode of measurement, you can enable a software emergency switch off of the PP- device (Option: **Enable Software Switch-OFF**). This emergency switch off should help to avoid a damage of your probes due to over- voltages and over- currents respectively. Therefore, insert the values (**Security conditions**) where the switch off is to be performed.

But note that a software switch off is not a guarantee because if the communication between the PC and the PP- device is lost, for instance due to a PC hang-up, the emergency switch off will fail.

**Maximum of current** (+) : determines the upper (anodic) current

Maximum of current (-)

Maximum potential (+)

Maximum potential (-)


: determines the lower (cathodic) current
: determines the upper (anodic) potential
: determines the lower (cathodic) potential

**FORCE RESET:** Clicking on this button, a software reset is performed. But keep in your mind that all measured data which are not saved will be lost.

Below the FORCE RESET button, you find a group box where you can select the maximum number of measured points per second. You can select one of the entries before you start a measurement. Please note that the selected value is more an index because the point density is essentially determined by the electrochemical parameters of each particular type of measurement (see below)

## 3.1 Dynamic measurements

The 'turnover' potentials / currents for the dynamic measurements are inserted within the table **Dynamic parameters** – on the left hand side of the figure above.

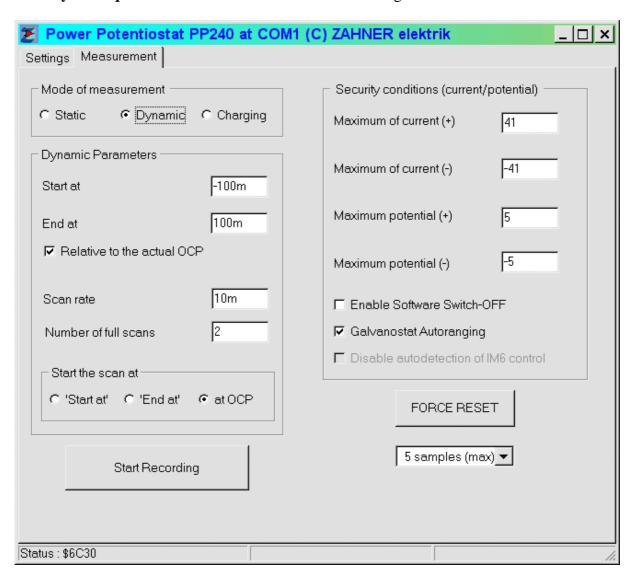Start at : defines the lower (= cathodic) turnover potential or current repectively.
 End at : defines the lower (= anodic) turnover potential or current repectively.

**Relative to the actual OCP:** defines whether the value above (**Start at** and **End at**) should be referred to the rest (or open circuit potential).

For instance (potentiodynamic mode): with **Start at** = -100 m, **End at** = +200 m and a measured rest potential of +250 mV results a lower turnover potential of +150 mV (absolute) and an upper turnover potential of +450 mV, provided the **Relative to the actual OCP** feature is selected. If not selected, the turnover potentials are set to -100 mV and +250 mV, independent of the rest potential.


**Start the scan at :** determines the point to start the dynamic measurement (scan).

**Scan rate**: determines the scan rate (in s<sup>-1</sup>). The smallest scan rate is 1 mV/s and 1 mA/s respectively. Provided in **Start the scan at** the option '**at OCP**' is selected, the sign of **Scan rate** determines whether the scan start to more cathodic values (negative scan rate) or to more anodic values (positive scan rate). In all other cases the sign is ignored and set appropriately.

**Number of scans:** determines the number of cycles to be performed. A cycle is defined as the 'range' between **Start at** and **End at.** The following table gives examples for different standard measuring techniques depending on **Start the scan at** and **Number of scans**:

Setting	Setting	Electrochemical
Start the scan at	Number of scans	technique
At OCP	2	Cyclovoltammetry (1 cycle)
At OCP	4	Cyclovoltammetry (2 cycles)
Start At	1	ramp (anodic to cathodic scan)
End At	1	ramp (cathodic to anodic scan)

Clicking on the button **Start Recording** will demand for a filename to save the measurement.

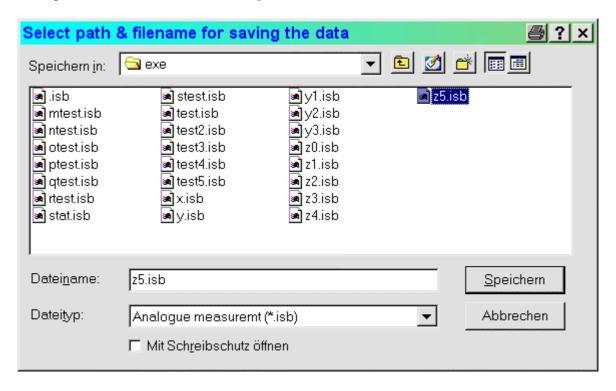After inserting a filename, the measurement will be started. There a two files, created when the measurement is finished: A binary file with the extension \*.isb (reserved for future use) and a file with the extension \*.txt which contains the measured values in ASCII format, ready to input in any calculation program.


#### Notes:

- 1.) Provided a measurement is running actually, you are asked to interrupt this measurement.
- 2.) If you don't define a filename completely, no measurement will be started.
- 3.) The procedure to start a dynamic measurement is as follows:
  - f. The PP- device is switched off.
  - g. You define a filename for saving the measurement.
  - h. The rest potential is measured.
  - i. The initial value will be set.
  - j. The PP- device is switched ON and the measurement will we started.

Therefore, switching ON or OFF is performed automatically.

# 3.2 'Charging' measurement

This type of measurement enables the galvanostatic charging and discharging of rechargeable systems like accumulators.

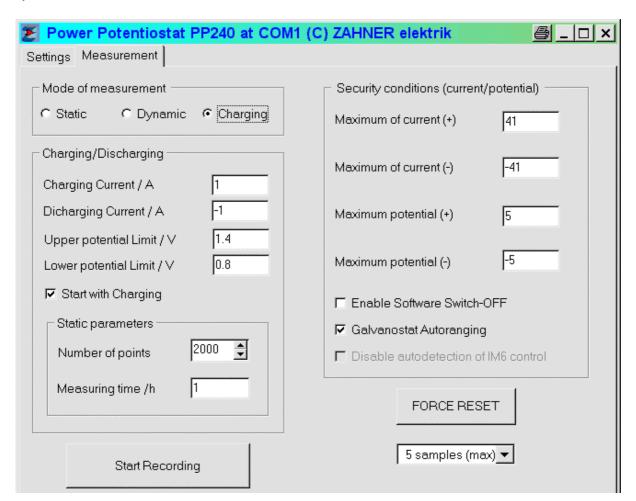
With the entries **Charging current** and **Discharging current** you determine the (static) currents you want to apply consecutively during the cycling of your system. Your system will be charged, until the **Upper potential limit** is reached. At this point, the **Discharging current** is applied until the **Lower potential limit** is reached, and so on.

Enabling the **Start with Charging**, the measurement will be started with the **Charging current**, otherwise the **Discharging current** is applied.

Additionally, you ca select the **Number of points** to be recorded as well as the overall **Measuring time**.

### 3.3 Static measurement

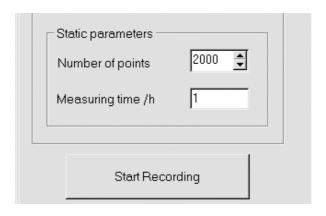The parameters to 'perform' a static measurement are set in the table **Static parameters** (right hand side of the figure above).

**Measuring time:** desired measuring time (in hours).

**Number of points :** (Approximate) Number of data points to be measured. The exact number of measured points will be determined from the time to be measured and the number of measured points per second.

The measurement will be started as described in the dynamic section except that the ON/OFF condition of the potentiostat will not be changed anyway.

# 4. Diagram

In this menu, a measurement is displayed graphically.


The diagram shown below depicts an actual recorded measurement. On the right side of the diagram you can select the diagram time to be displayed:

5.) current vs time: Plot of the current as a function of time6.) potential vs time: Plot of the potential as a function of time

7.) **current vs potential**: Plot of the current as a function of the potential 8.) **potential vs current**: Plot of the potential as a function of the current


Pressing the secondary mouse button (usually the right one) in the diagram will offer the possibility to export the graphics immediately. For instance, you can copy the actual graphics into the windows clipboard or export it as an EMF- graphics file.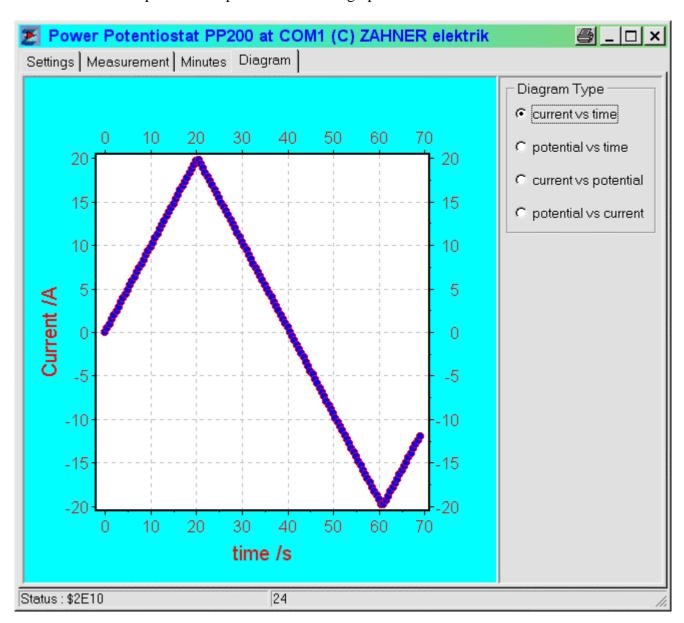

### 5. Minutes

Starting a measurement, the most important parameters are noted in the text- field **Minutes** . You can add own comment if you like as long as the measurement is running. You can 'save' the minutes simply by a 'copy and paste operation' into any text document you like.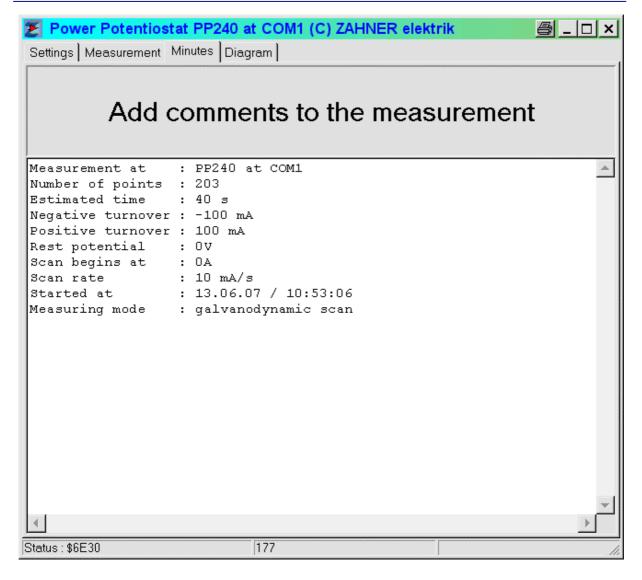

Note: As long as you have not started a new measurement, the minutes will be updated regardless whether the measurement has been stopped.